

# CMEB-REF-ST-12-G-NSF

### **Product Description**

These cutting-edge pharmacy refrigerators are certified in accordance with the NSF/ANSI 456 Standard for Vaccine Storage. With this certification, units protect pharmaceuticals at optimal temperatures, preventing waste and allowing for peak delivery. Our Standard line provides multi-functional features in a cost-effective design.

These glass door refrigerators utilize microprocessor controllers and feature temperature alarms, remote alarm contacts, and probe access ports with included probes. Units run on natural, hydrocarbon refrigerant for environmental health and energy efficiency.

#### **General Description and Application** Description Single Glass Door Pharmacy/Vaccine Upright Refrigerator Operational environment Indoor use only, +18°C to +26°C (+65°F to +78°F), <70% RH Storage capacity 12 cu. ft. gross volume Door One swing glass door, self-closing, right hinged, non-reversible, magnetic sealed gasket, keyed Four shelves (three adjustable/one fixed) with guard rail on back Shelves 3 1/2" Swivel Casters(two locking) Mounting Shielded, switched LED lighting, full coverage, balanced spectrum Interior lighting Forced Air technology, patent pending Airflow management Rear wall port (3/4") dia. External probe access Cabinet is foamed-in-place with EPA compliant high density urethane foam Insulation **Exterior materials** White powder coated steel Access control Pyxis®, Omnicell® and AcuDose RX® compatible One (1) year parts and labor warranty, excluding display probe calibration General warranty Compressor warranty Five (5) years compressor warranty 224 lbs. **Product Weight** Shipping Weight 264 lbs. Rated Amperage NEMA 5-15 plug, 8 to 10 ft typical, conforms to UL471 requirements, Vaccine Storage power Power Plug/Power Cord 110-120V AC: 15 A (minimum) Facility Electrical Requirement Agency Listing and Certification Compliant with the temperature performance requirements as defined in the NSF 456 Standard for Vaccine Storage for all testing scenarios. UL, C-UL, ETL, C-ETL listed and certified to UL471 standard, hydrocarbon refrigerant safety. Pharmacy refrigerator/freezer toolkit and temperature logs **Included Accessories**

## Refrigeration System

Compressor

Refrigerant

Condenser

Evaporator

Evaporator

Hermetic, high performance

EPA SNAP compliant, R290, propane

Fin and tube design, high efficiency fan

Fin and tube design, high efficiency fan

Cycle optimized, zero energy

Performance	
Uniformity <sup>1</sup> (Cabinet air)	+/- 0.7°C
Stability <sup>2</sup> (Cabinet air)	+/- 1.3°C
Maximum temperature variation	+/-1.1°C
(Cabinet air)	
Temperature rise after 8 sec door	Temperature did not exceed 5.4°C at any probe for all required NSF/ANSI 456 testing
openings	protocols <sup>3</sup>
Recovery after 3 min door opening	All probes recover to under 8°C within 2.5 min.
Energy consumption	0.61 KWh/day⁴
Average heat rejection	1.20 KWh/day (315 BTU/h)⁴
Noise pressure level (dBA)	48 or less installed
Pull down time to 4°C nominal operating	30 min
temn	

::-::p		
Controller, Configuration, Alarms	and Monitoring	
Controller technology	Parametric, microprocessor, LED display with 0.1°C resolution	
Temperature setpoint range	1°C to 10°C (Controller settings must remain unaltered to ensure thermal performance	
	compliant with NSF/ANSI 456 Standard for Vaccine Storage requirements)	
Calibration	Calibrated using a NIST traceable device, certificate included	
External alarm connection	State switching remote alarm contacts	
	Visual and audible indicators	
Alarms	High / Low temperature, compliant with alarm requirements defined in the NSF/ANSI 456	
	Standard for Vaccine Storage	
Simulator ballast	Glass bead thermal media	

Performance data acquired at 22°C ambient, using NSF/ANSI 456 compliant validation ballast probes, empty chamber, during stabilized steady state operation and a DAQ sampling rate of one measurement every 10 seconds

- 1 Uniformity is defined as the maximum variance in temperature across all probes at any point in time over the testing period
- 2 Stability is defined as the maximum variance in temperature experienced by any single probe over the testing period
- 3 Temperature performance for all loaded and unloaded door opening protocols, all alarm, controller and probe requirements as defined in the NSF/ANSI 456 standard for vaccine storage
- 4 Data per Energy Star test results or equivalent testing and calculation. Heat rejection based on daily averages, not continuous operation. Performance exceeds Energy Star requirements.

#### **Product Data Sheet**

Upright 12 cu. ft. Glass Door Refrigerator, High Performance - Certified to NSF/ANSI 456 Standard for Vaccine Storage


#### Certifications








\*-one or more of these certifications may apply to this unit.

Temperature Probes								
Probe	Ave Min		Max					
1	3.5	2.7	4.7					
2	3.8	3.4	4.5					
3	4.0	3.7	4.5					
4	3.7	3.1	4.7					
5	3.8	3.4	4.5					
6	3.9	3.4	4.7					
7	3.8	3.4	4.6					
8	4.3	3.9	5.0					
9	3.5	2.7	4.8					
10	3.8	3.2	4.7					
11	3.6	3.1	4.6					
12	3.6	3.1	4.4					
13	3.7	3.4	4.3					
14	4.2	3.8	4.9					
15	3.4	2.8	4.5					

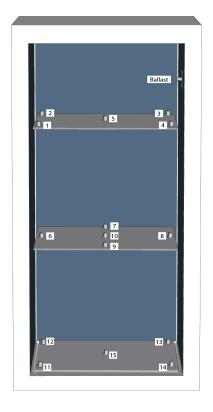

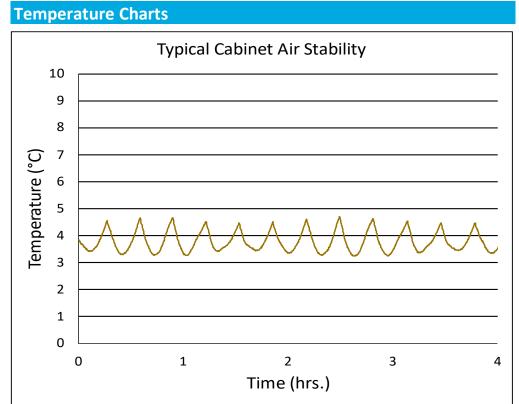

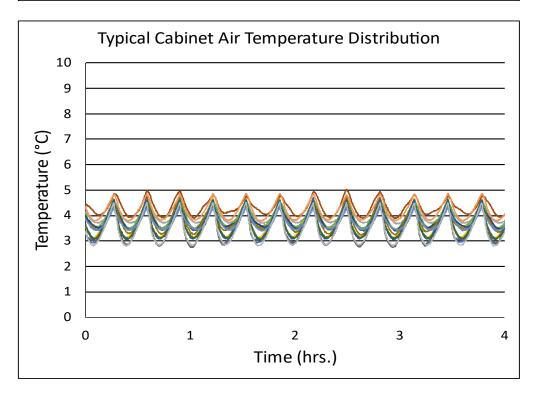

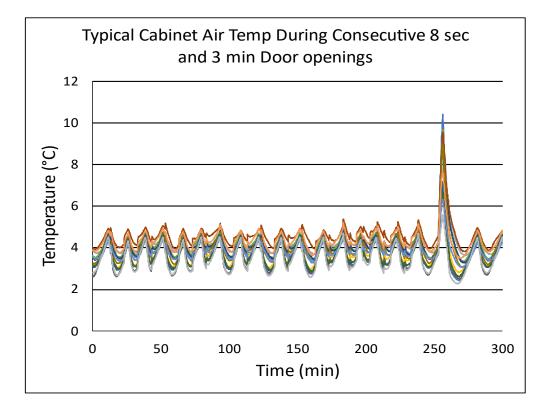



### **Product Data Sheet**

Upright 12 cu. ft. Glass Door Refrigerator, High Performance - Certified to NSF/ANSI 456 Standard for


Vaccine Storage


### **Images**





Dimensions								
	Width	Depth	Height	Door Swing	Total open Depth			
Exterior	25"	29"	65 3/4"	23 1/4"	52 1/4"			
Interior	20 3/8"	19 1/4"	36 1/8"					

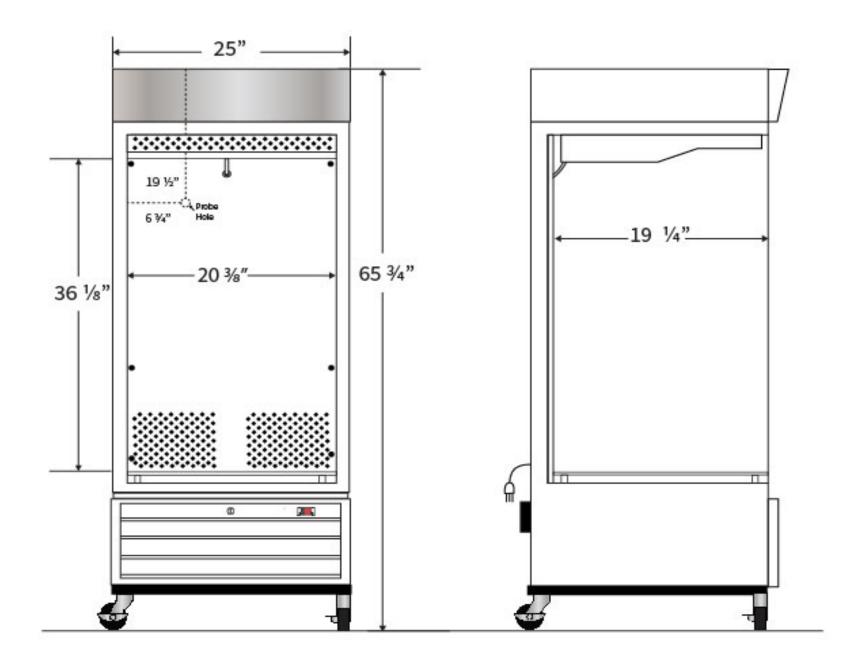

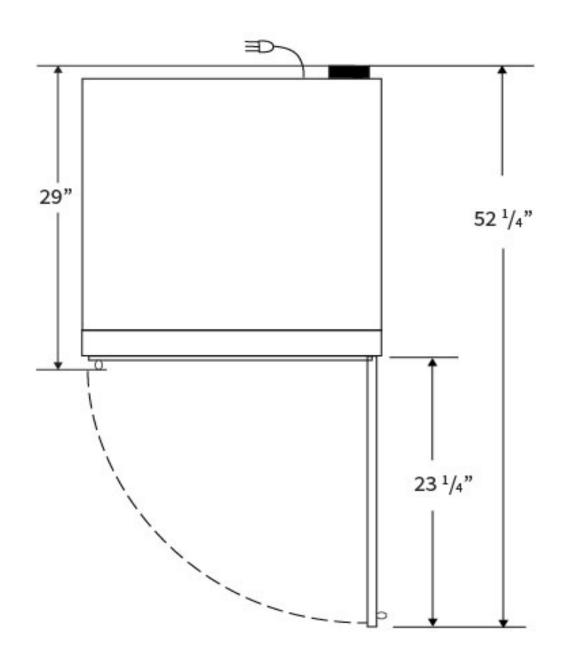

Contact
Customer Serv

800.338.2372

customerservice@cmecorp.com